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3-jaw parallel self-centering pneumatic gripper (series T)

- Modular with Gimapick system.
- Double acting.
- Air feeding possible directly from the fixing plate.
- Possibility to mount it on front with through screws.
- High gripping force with low weight.
- Optional magnetic sensors.

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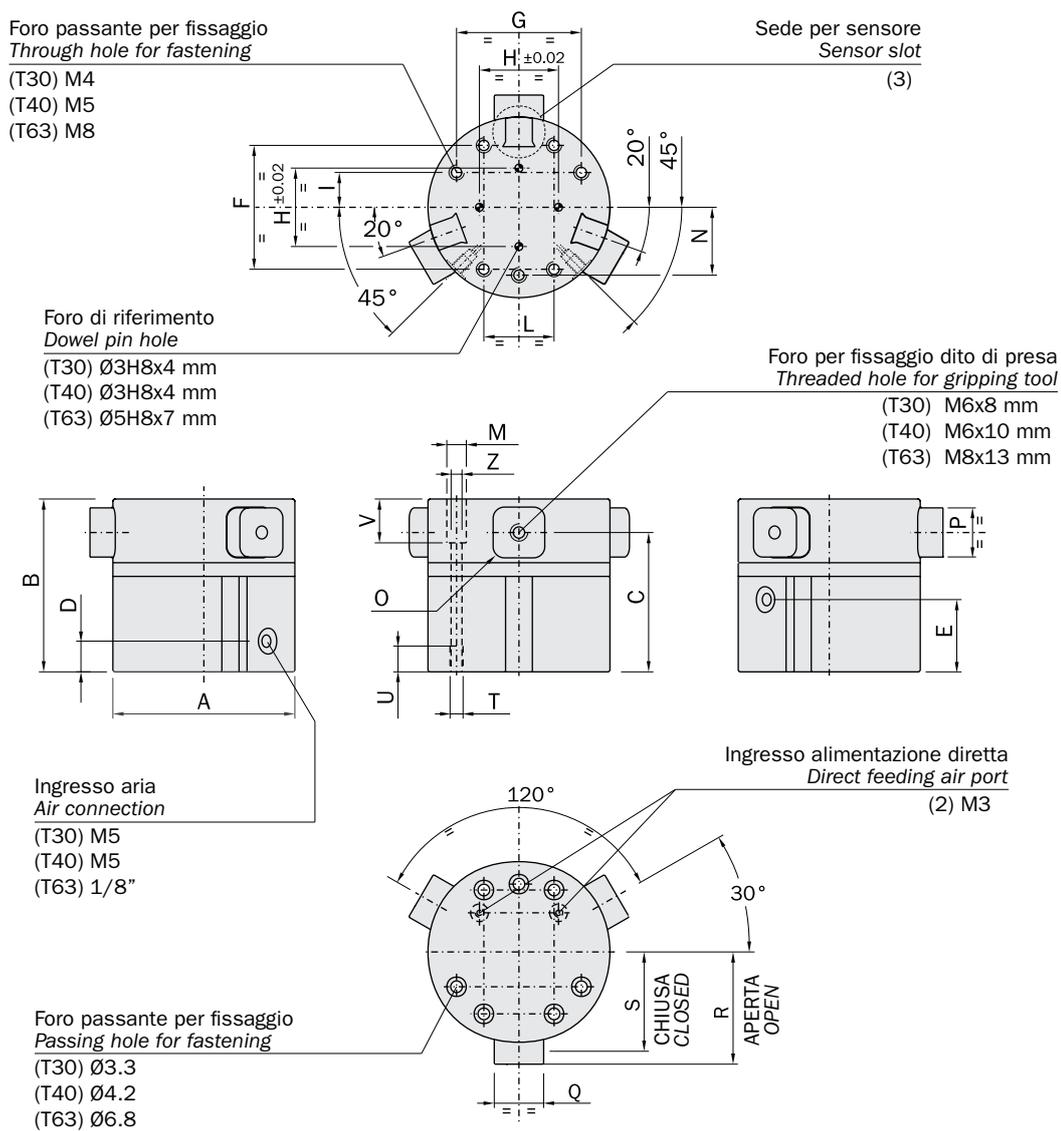
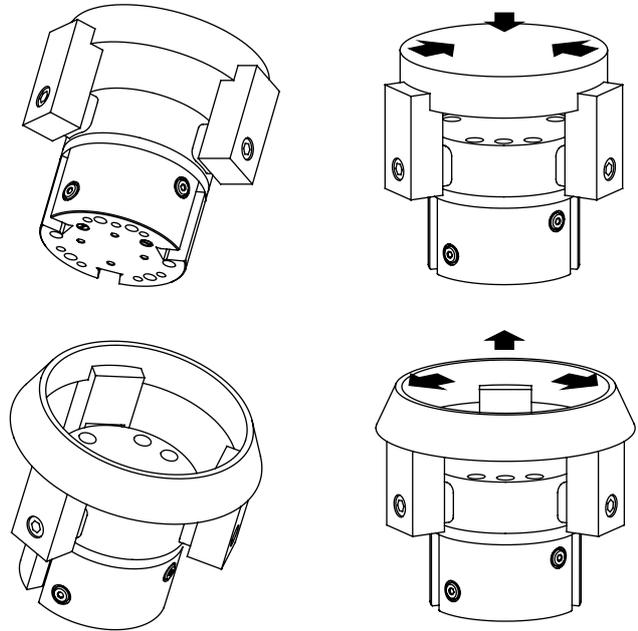
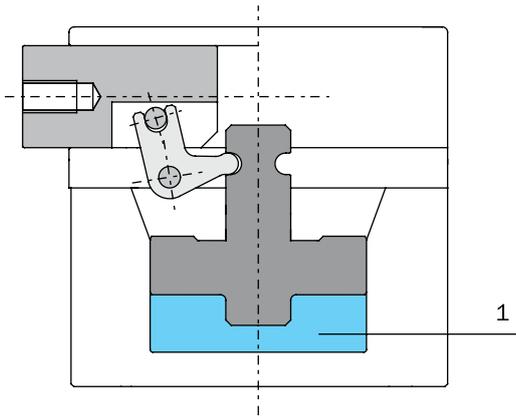
Gripping

The three jaw gripper allows the correct gripping of pieces with different diameter keeping the center in the same position. The jaws are linked to three levers moved by one piston, having by this, self-centering action.

The gripper is double-acting, so internal or external gripping is possible.

Camera in pressione
Pressurized chamber

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

Application examples

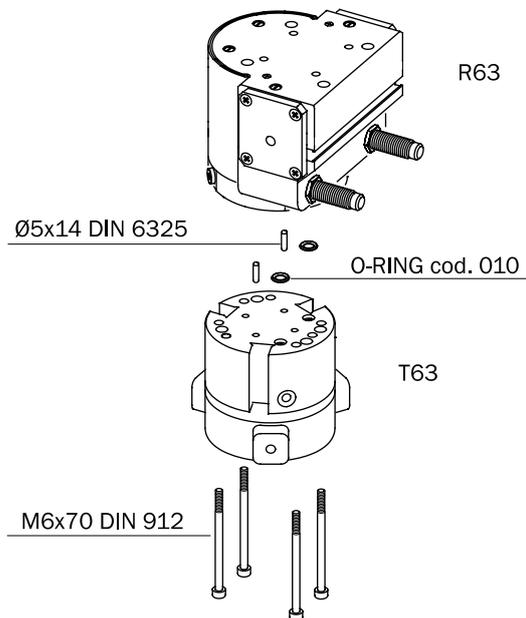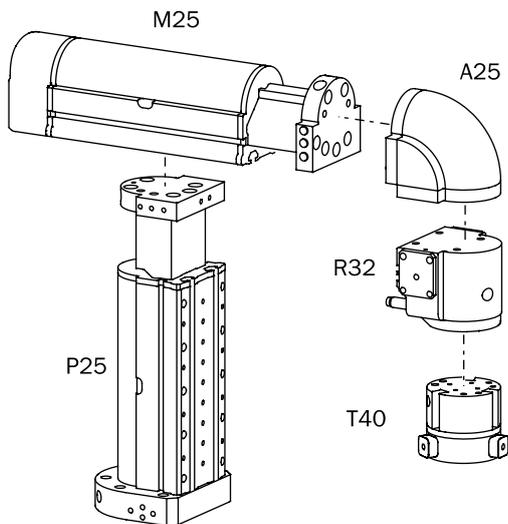These grippers can be easily mounted with Gimapick elements.

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


Gripping force

The graphs show the gripping force on each jaw, as a function of the operating pressure and the gripping tool length (Z).

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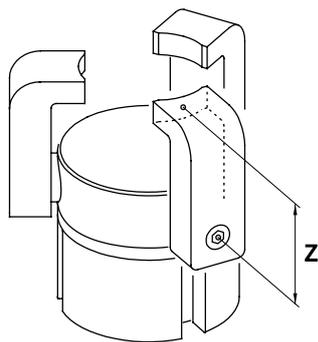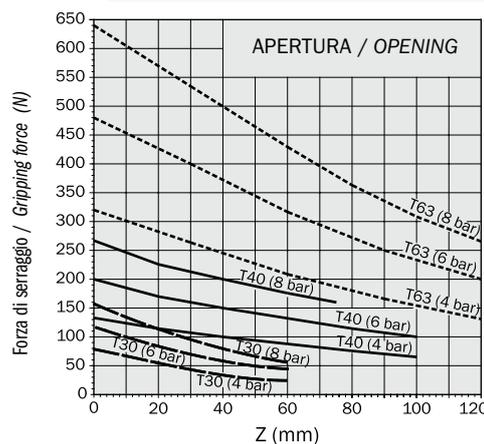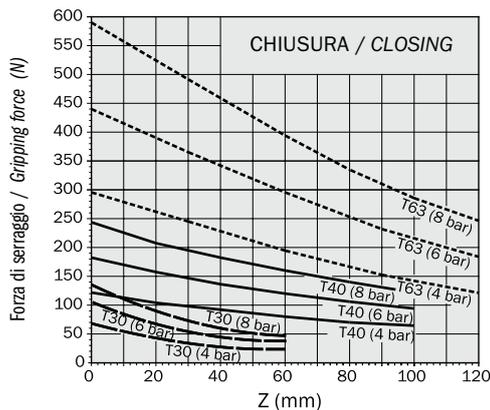The force shown in these graphs refers to one jaw. The total force is triple.

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
Safety loads

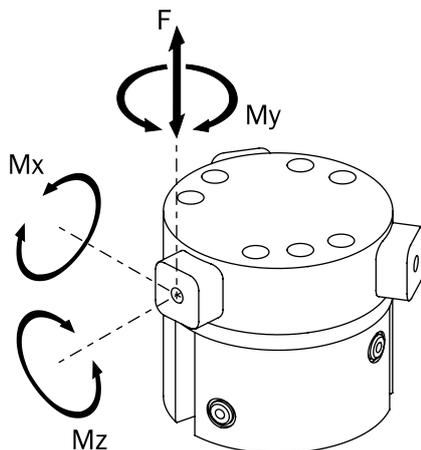Check the table for maximum permitted loads. Excessive forces or torques can damage the gripper, cause functioning troubles and endanger the safety of the operator. F s, Mx s, My s, Mz s, are maximum permitted static loads. Static means motionless jaws. F d, Mx d, My d, Mz d, are maximum permitted dynamic loads. Dynamic means running jaws. The following table shows the specified maximum loads (m) on each gripping tool as a function of closing or opening time. Use flow controllers (not supplied) to get the proper speed.


	T30	T40	T63
F s	100 N	300 N	500 N
Mx s	4 Nm	12 Nm	32 Nm
My s	4 Nm	12 Nm	32 Nm
Mz s	4 Nm	12 Nm	32 Nm
F d	1 N	3 N	8 N
Mx d	3 Ncm	8 Ncm	30 Ncm
My d	3 Ncm	8 Ncm	30 Ncm
Mz d	3 Ncm	8 Ncm	30 Ncm
m 0.2 s	150 g	300 g	700 g
m 0.1 s	100 g	200 g	500 g
m (*)	80 g	150 g	400 g

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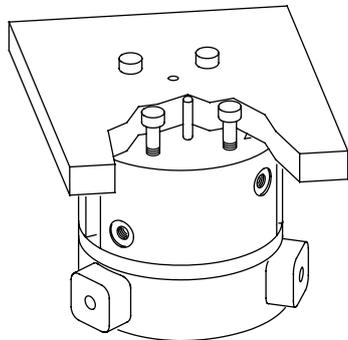
Gripper fastening

The gripper can be fastened to a static or moving part.

When on a moving part, you must pay attention to the forces created by inertia on the gripper and its load.

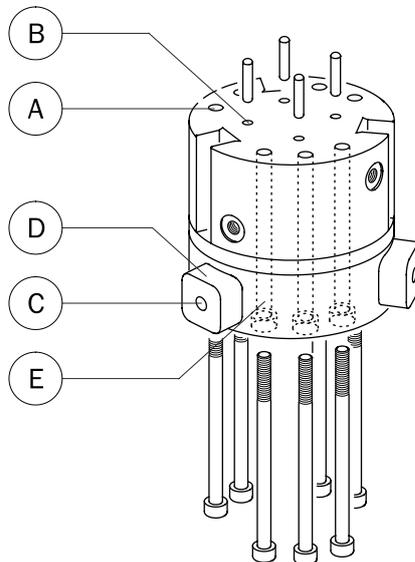
To fasten the gripper use at least four screws and two dowel pins.

	T30	T40	T63
A	M4x8 mm	M5x10 mm	M8x16 mm
B	Ø3H8x4 mm	Ø3H8x4 mm	Ø5H8x7 mm
E	Ø3.3 mm	Ø4.2 mm	Ø6.8 mm

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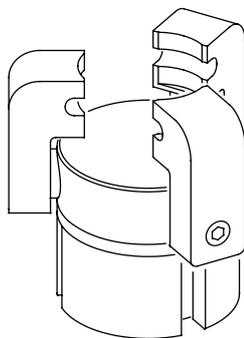
Gripping tool fastening

The gripping tools must be as short and light as possible.

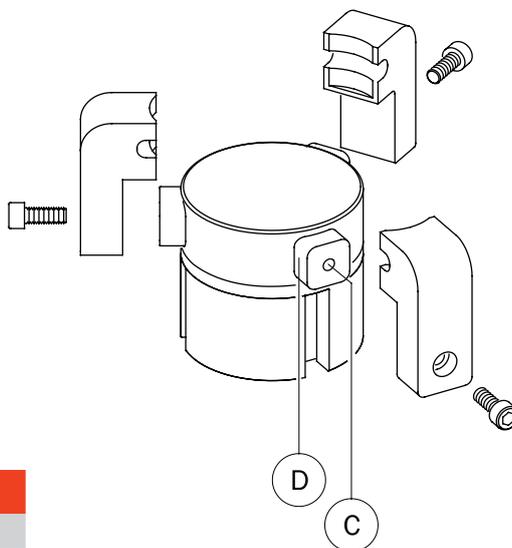Fit them to the jaws by centering the square calibrated profile (D) and locking with a screw through the threaded middle hole (C).


	T30	T40	T63
C	M6x8 mm	M6x10 mm	M8x13 mm
D	14 ^{-0.05} x 14 ^{-0.05} mm	19 ^{-0.05} x 19 ^{-0.05} mm	26 ^{-0.05} x 26 ^{-0.05} mm

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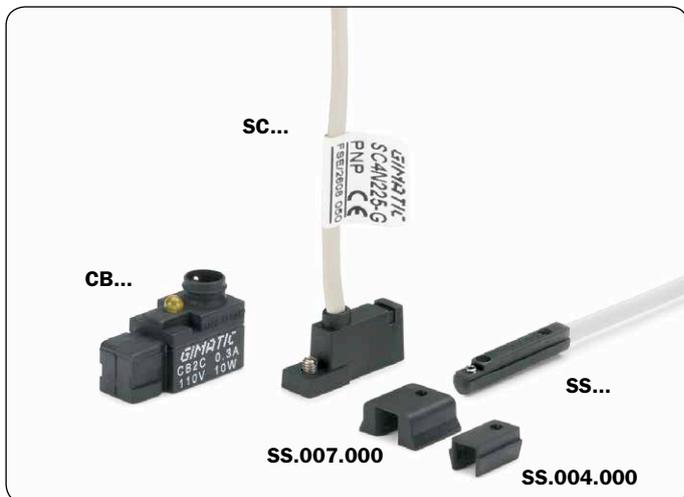
Sensors

The operating position is detected by proximity magnetic sensors (optional) through a magnet placed on the piston. Therefore, avoid using the gripper in the vicinity of intense magnetic fields or near a large mass of ferromagnetic material as this may cause detection errors.

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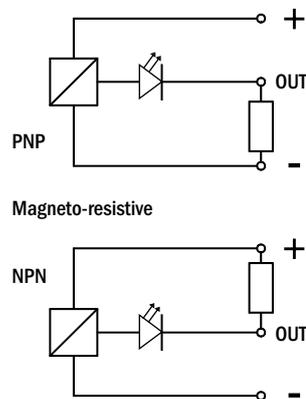
(1)
By the adapter (SS.007.000) provided with the pack K-SENS.

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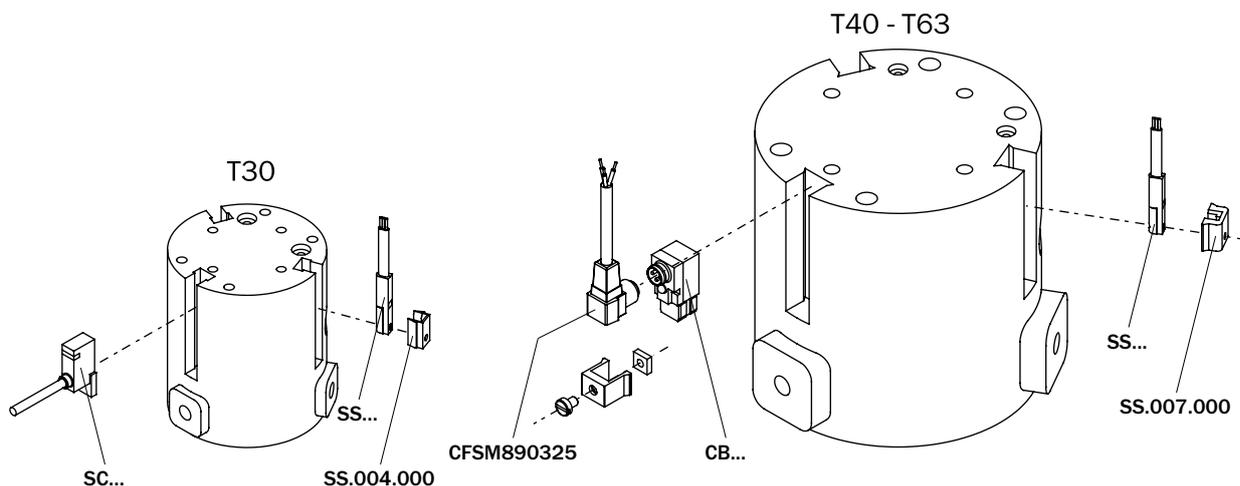(2)
By the adapter (SS.004.000) provided with the pack K-SENS.

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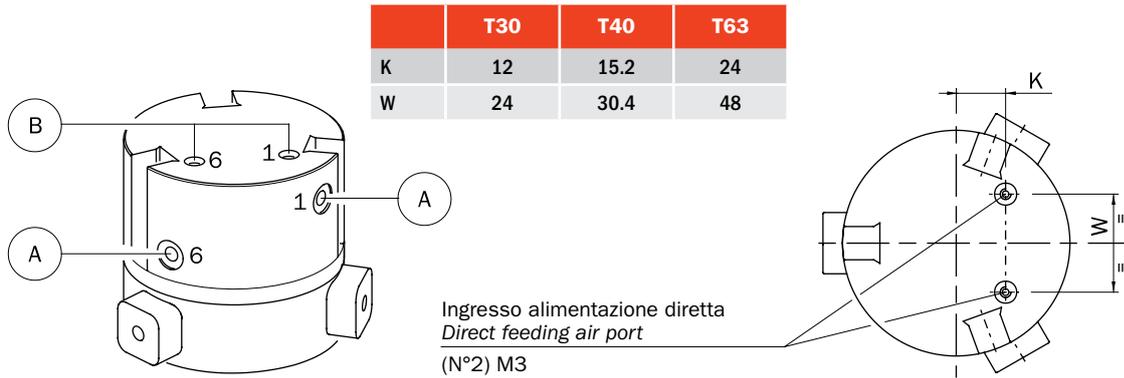
Compressed air feeding

The compressed air feeding is accomplished on the lateral air ports (A) with fittings and hoses (not supplied), or directly on the bottom air ports (B).

Compressed air in 1: gripper opening.
Compressed air in 6: gripper closing.

The compressed air, must be filtered from 5 to 40 µm.
Maintain the medium selected at the start, lubricated or not, for the complete service life of the gripper.

The pneumatic circuit must be pressurized progressively, to avoid uncontrolled movements.

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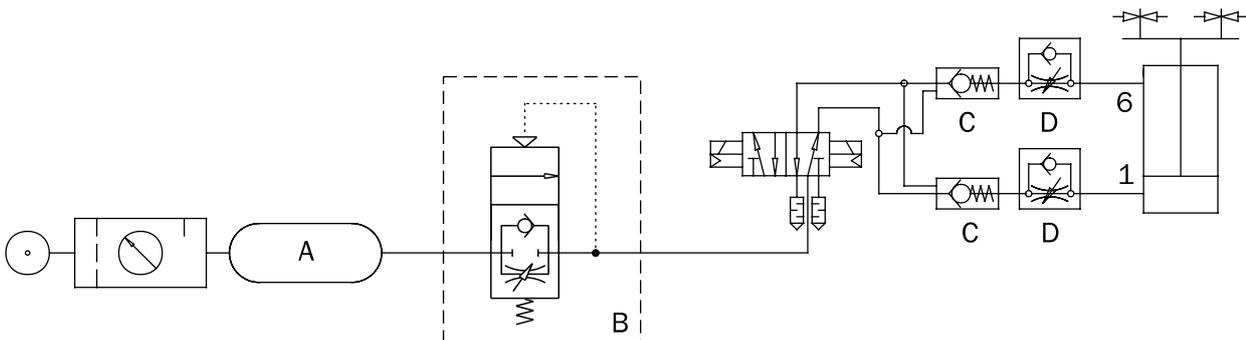
Pneumatic circuit

Possible problems on a compressed air circuit:

- 1- Pressure variation.
- 2- Pressurizing with empty cylinder.
- 3- Sudden pressure black-out.
- 4- Excessive speed of the jaws.

Possible solutions:

- 1- Compressed air storage (A).
- 2- Start-up valve (B).
- 3- Safety valve (C).
- 4- Flow controller (D).

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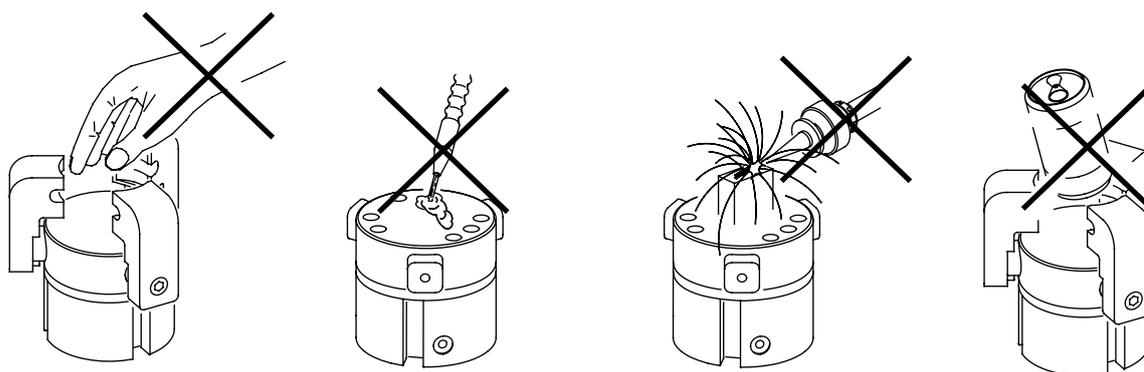
Caution

Never let the gripper come into contact with corrosive substances, soldering splashes or abrasive powders as they may damage the gripper.

Never let non-authorized persons or objects stand within the operating range of the gripper.

Never operate the gripper if the machine on which it is fitted does not comply with safety laws and standards of your country.

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Maintenance

Grease the gripper after 10 million cycles with:

- Molykote DX (metal on metal).
- Molykote PG75 (gaskets).

The figure below shows the jaw backlash.